

LABVIEW INSTALLATION INSTRUCTION

(AVAILABLE FOR WINDOWS ONLY)

[For MAC and Linux, kindly contact 5997]

STEP 1:

LICENSE REQUEST

- Before Install kindly copy the below link and paste it on the browser and fill the form to use this software. <http://web.iitm.ac.in/licenserequest/>
- To download the software, to login <https://hpce.iitm.ac.in/> using your LDAP credentials.
- After the login, on the right side below the **Quick links** → click **commercial software** to download the software.
- Download 2017LV-64WinEng.exe
- For remaining version contact HPCE team 5997
- Activation link for Labview https://www.ni.com/nlg-activate/jsp/customer_activate_details.jsp?p_serial_number=M21X71242

INSTALLATION

FOR WINDOWS:


STEP 2: Use **WINRAR** to extract the file

STEP 3: Run the setup as run as administrator

STEP 4: Click “Next”.



STEP 5: Click “Next”

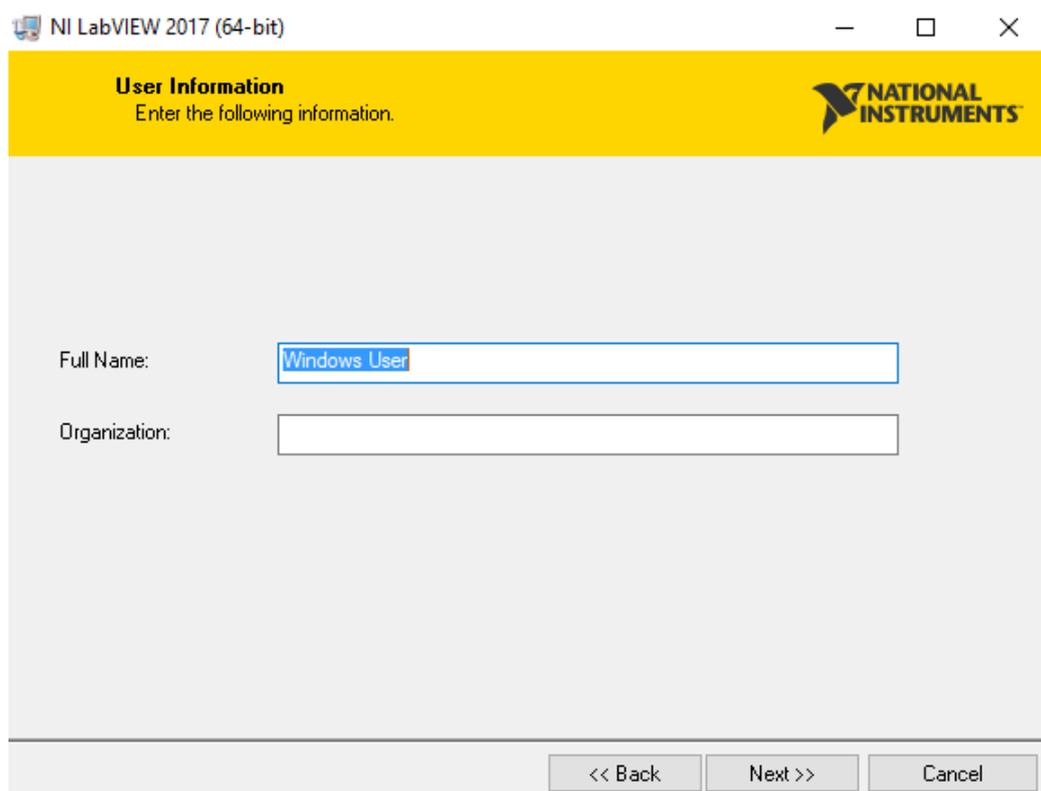

STEP 6: Click “Next”

NI LabVIEW 2017 (64-bit)

Serial Numbers
Enter Serial Numbers for the following Products

NATIONAL INSTRUMENTS

Serial Number:

LabVIEW 2017 (Base/Full/Professional) **

Application Builder for LabVIEW 2017 - leave blank if activating LabVIEW Professional. **

Report Generation Toolkit for LabVIEW 2017 - leave blank if activating LabVIEW Professional. **

**Leave serial number field empty to use product in evaluation mode.

<< Back Next >> Cancel

STEP 7: Click “Next”

NI LabVIEW 2017 (64-bit)

Destination Directory
Select the installation directories.

NATIONAL INSTRUMENTS

Select the folder to install NI Software


C:\Program Files\National Instruments\ Browse...

Select the folder to install NI LabVIEW 2017


C:\Program Files\National Instruments\LabVIEW 2017\ Browse...

<< Back Next >> Cancel


STEP 8: Click “Next”

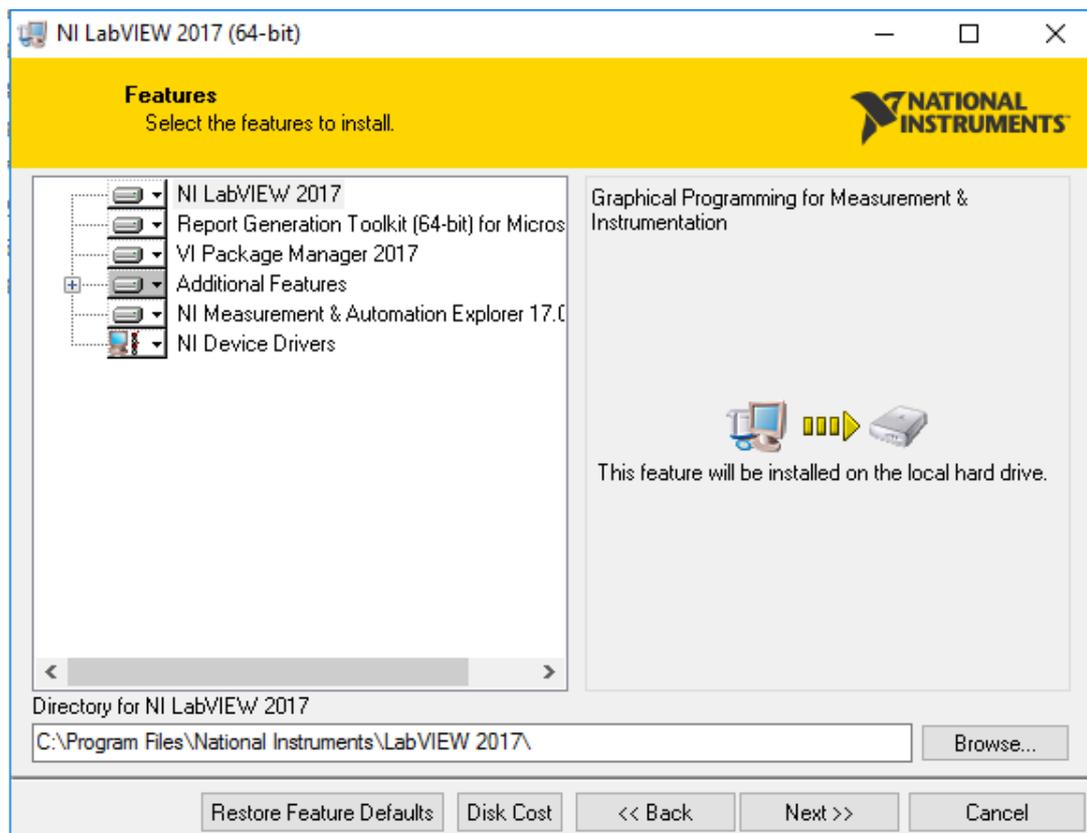


STEP 9: Click “Next”

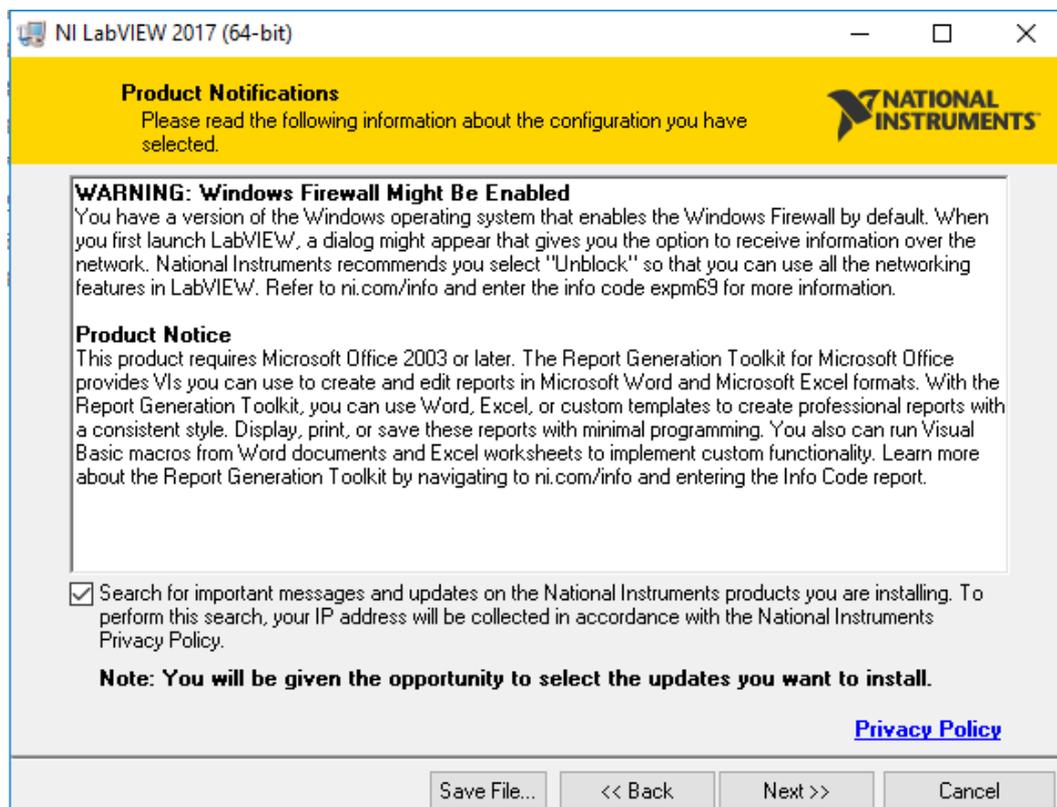


STEP 10: Click “Next”

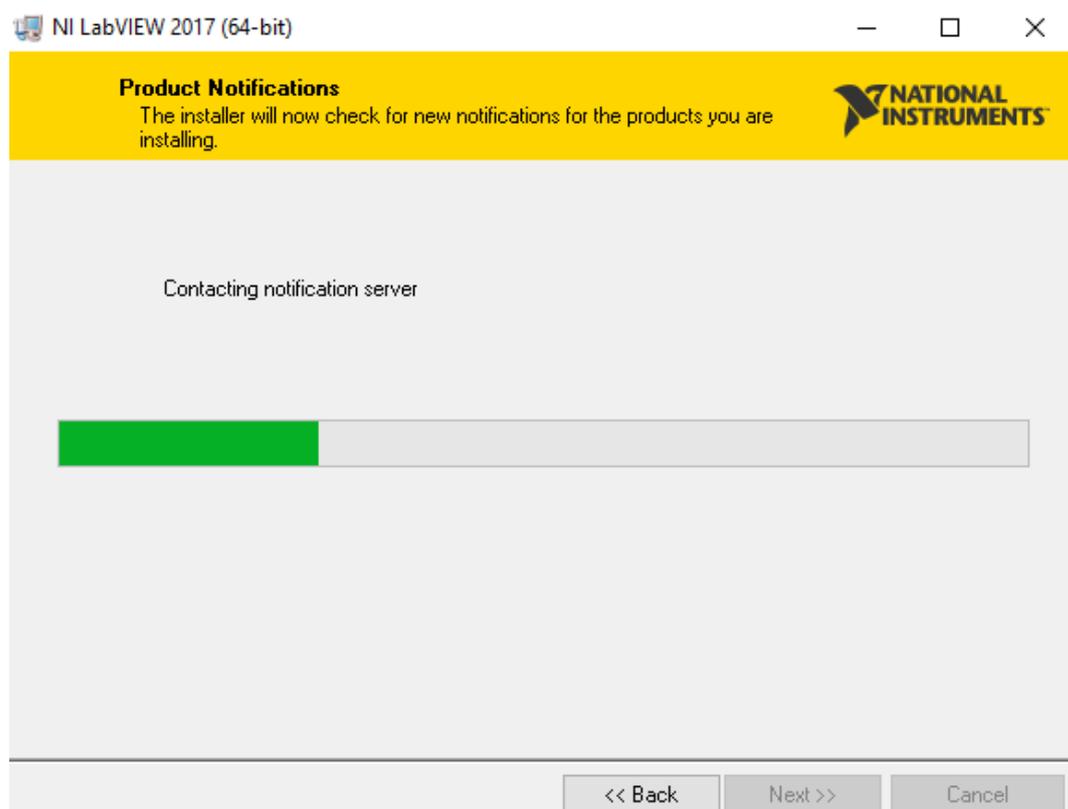


STEP 11: Click “Next”

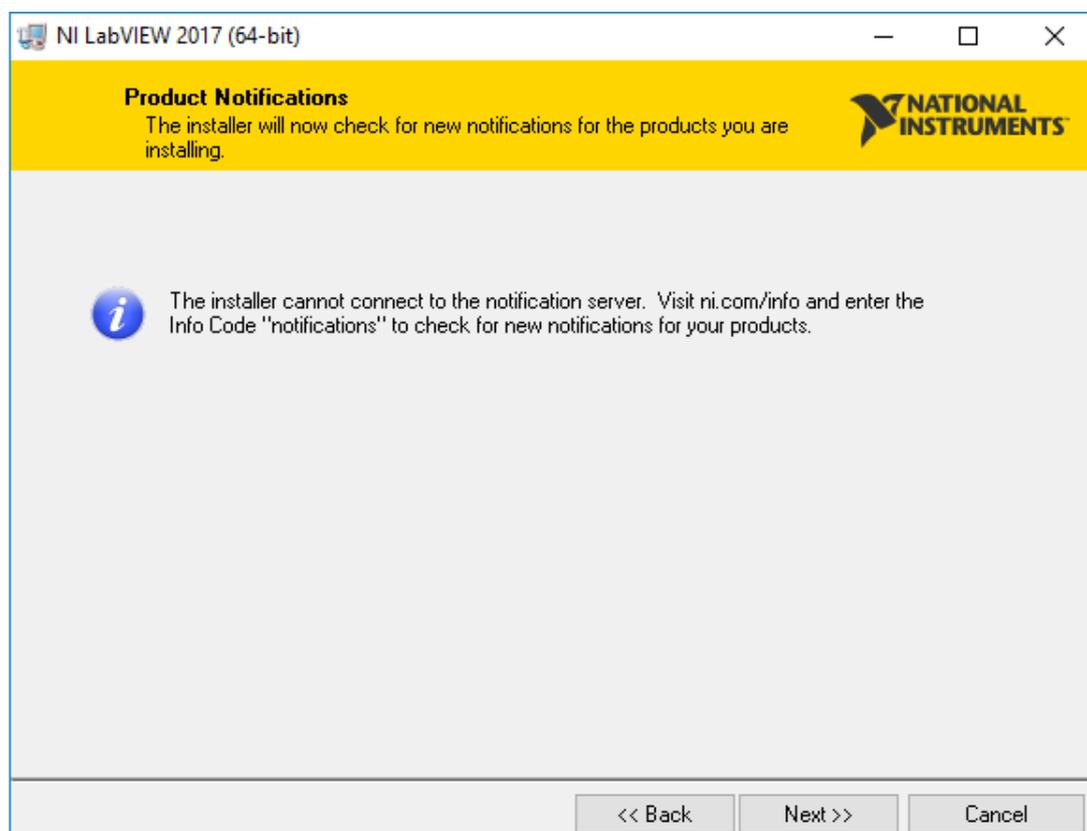


STEP 12: Click “I accept” & Click “Next”

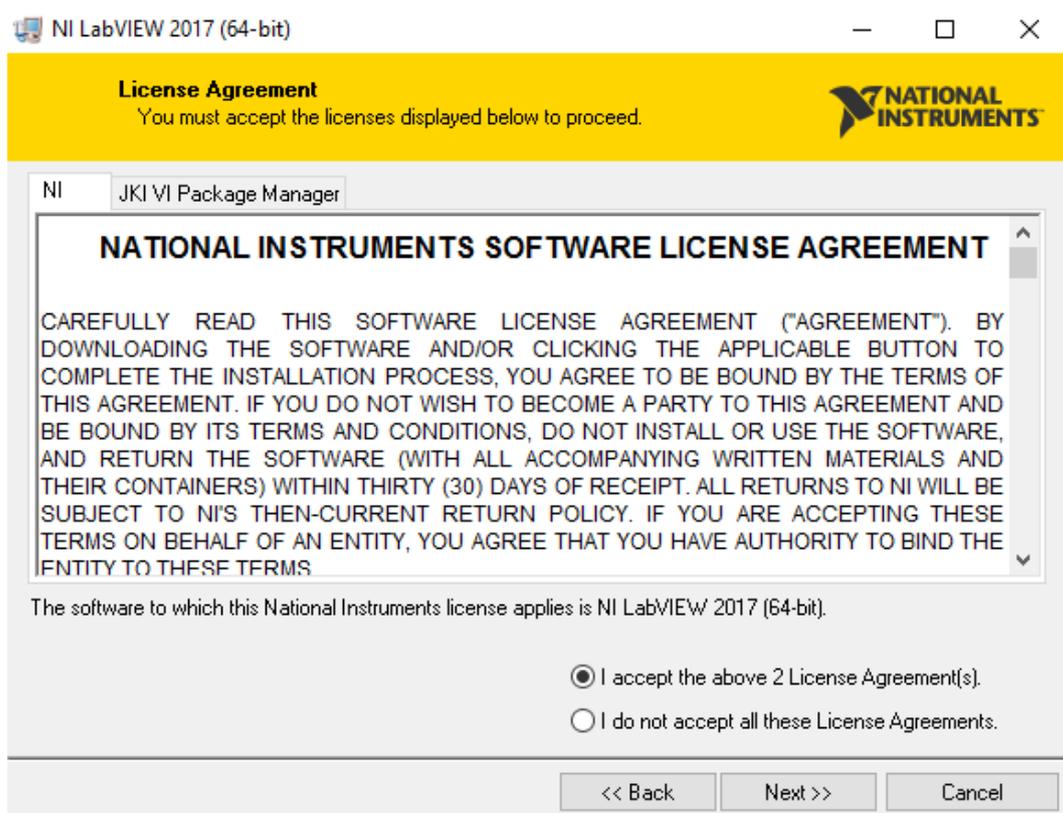


STEP 13: Click “I accept” & Click “Next”

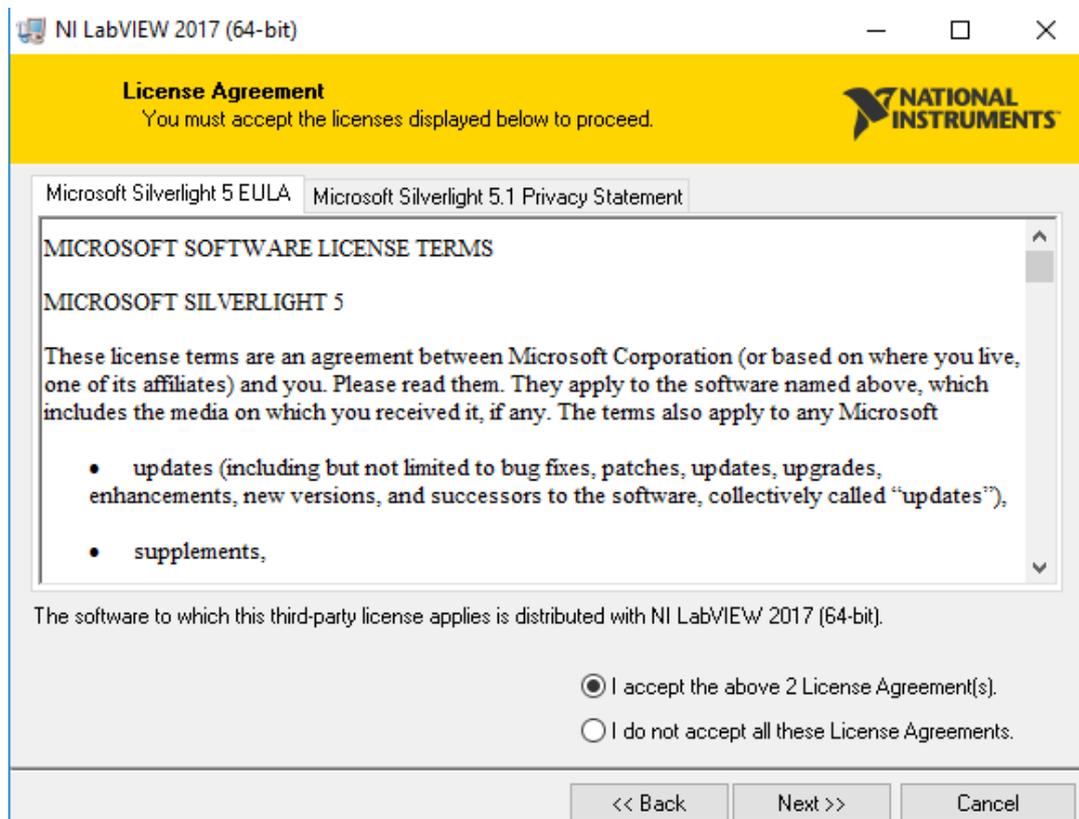


STEP 14: Click “Next”

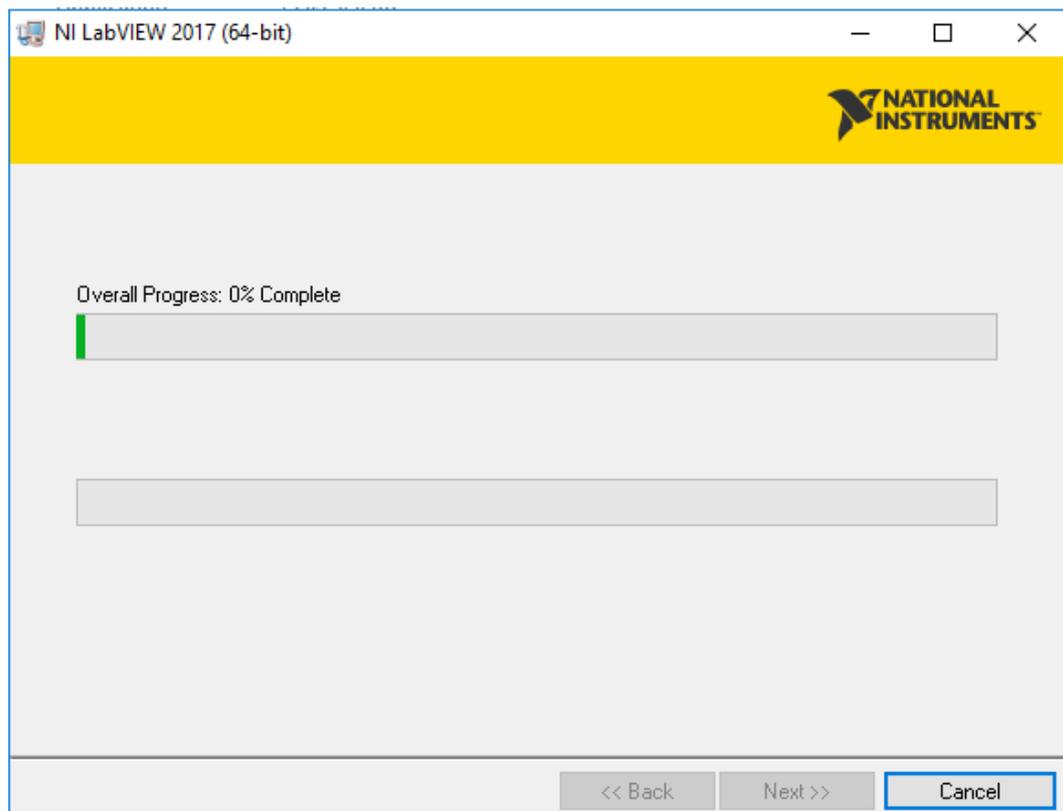

STEP 15: Click “Decline Support”

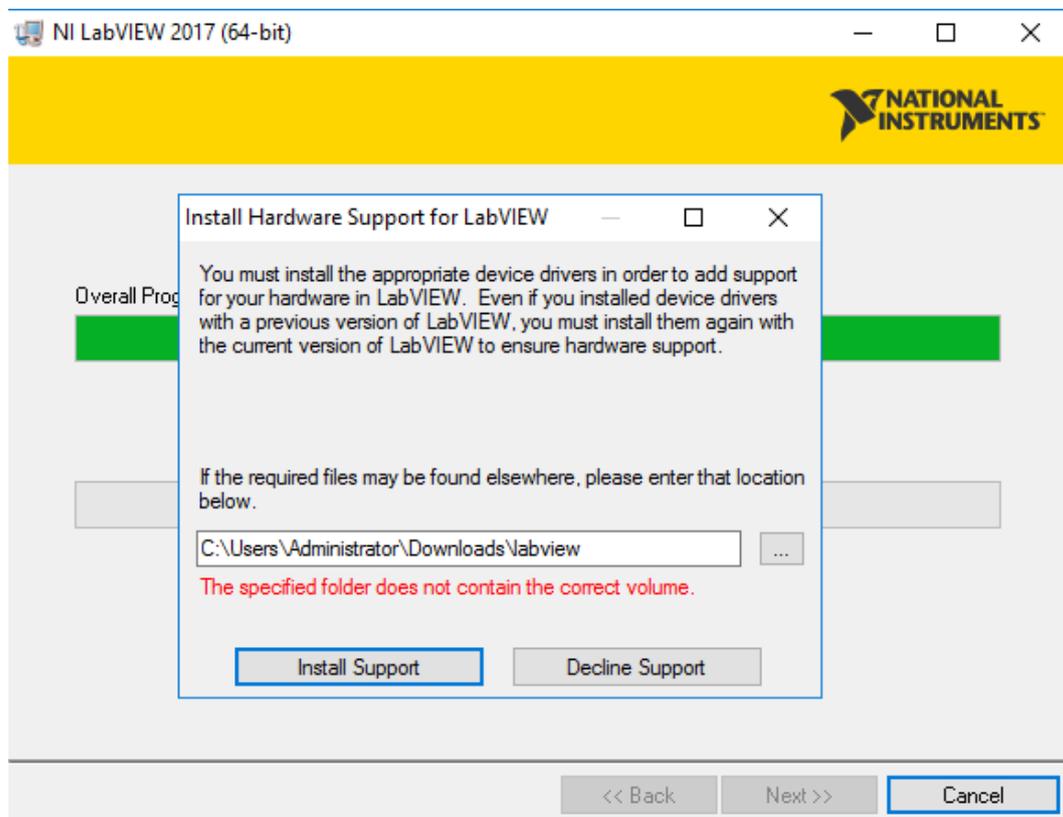


STEP 16: Click “Next”

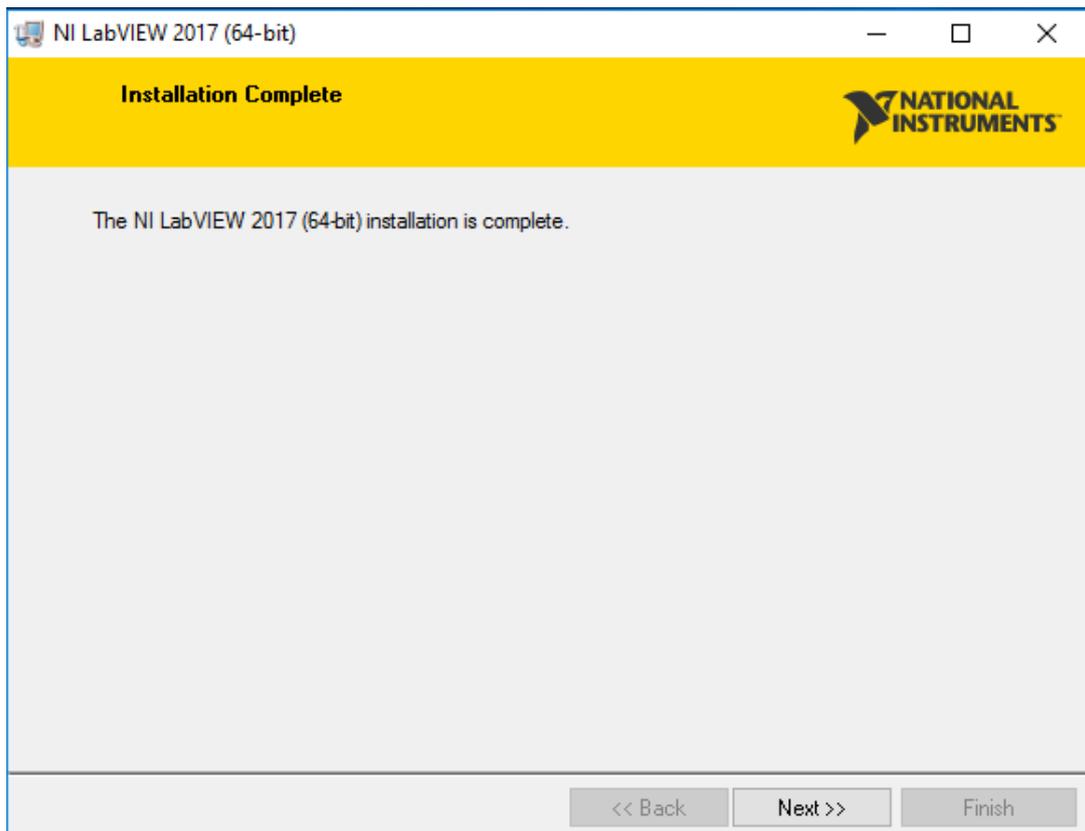


STEP 17: “Open LabView”


STEP 18: Click “Activate”

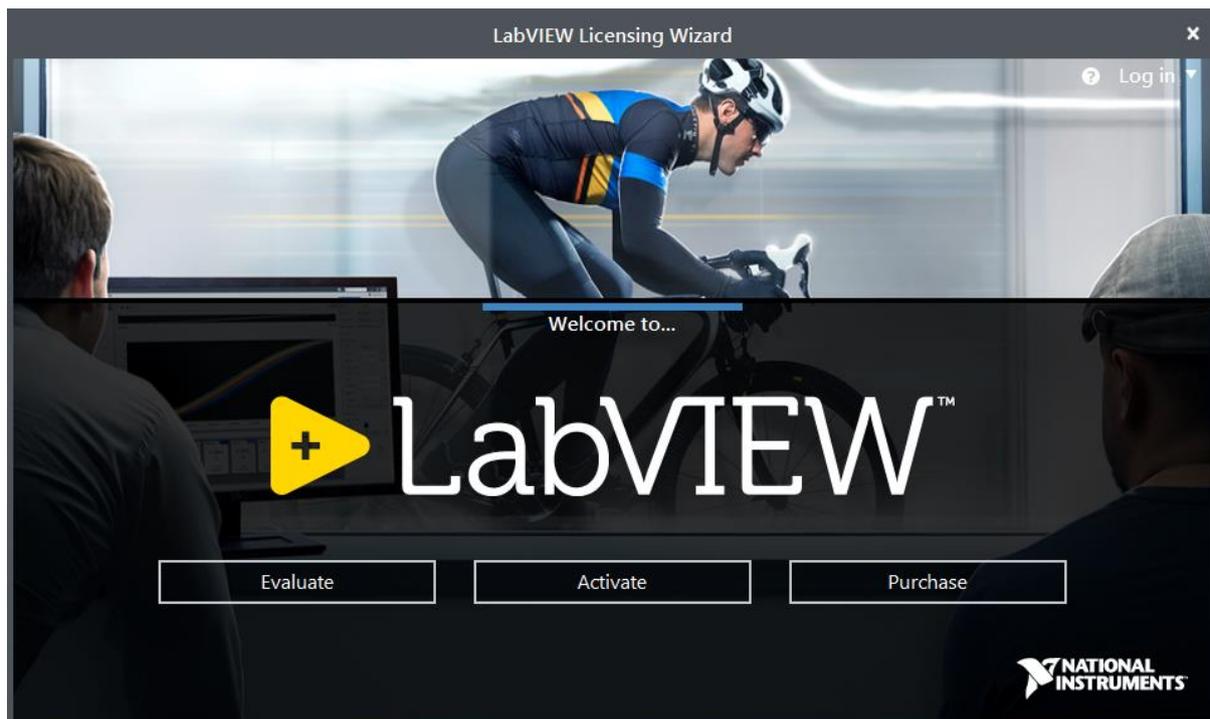

STEP 19: Click “X MARK”

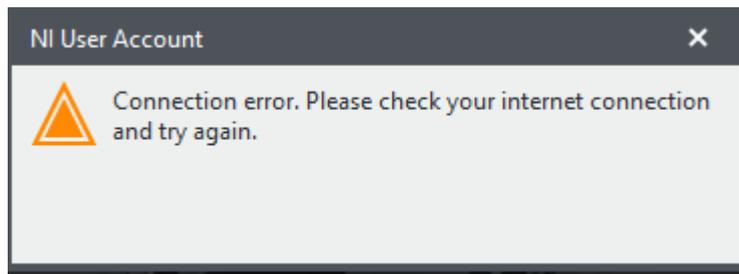

STEP 20: Copy the Activation Codes from the Activation link and paste here.

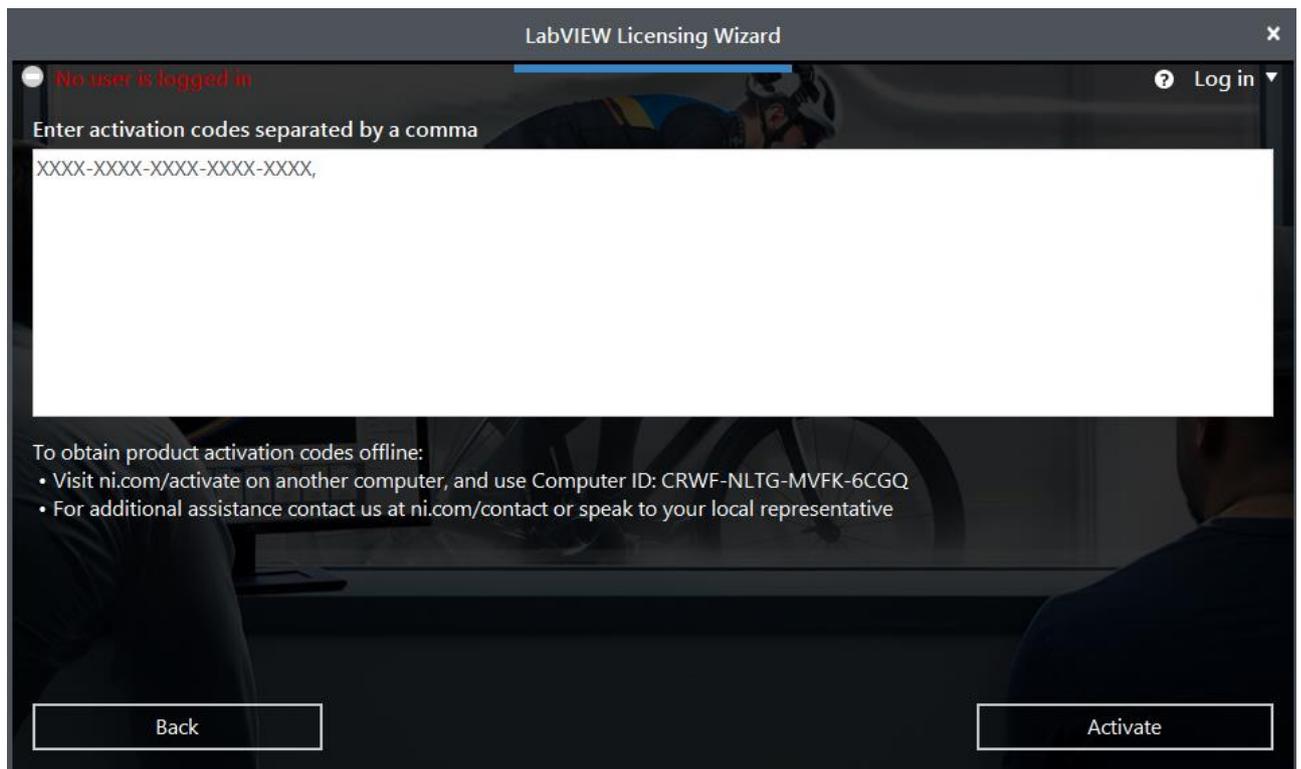

Now, the installation will be getting completed.